

Follow these steps to properly use your Full Kit 2018

1. **Lenses** – the Blips Full Kit 2018 includes four lenses: Macro Plus (yellow), Macro (green), Micro (blue) and Ultra (red), which respectively provide a magnification of 5x, 10x, 20x and 33x. Due to their high magnification rate, Micro and Ultra lenses need the supporting stage to provide their best performance.
2. **Lens on smartphone** – Stick the lens, well centered onto the smartphone camera (on the main camera for dual-cam phones). Micro and Ultra lenses: place the smartphone, with your Blips well aligned onto the camera, on the stage.
3. **Supporting stage preparation** - Remove the protective film from the stage parts. Mount the two equal rectangular plates when using the Micro lens on the smartphone. Slot the thicker rectangular plate (4mm-thick) with a little inclination to support the glass slide, when using the Ultra lens. Use the rubber bands around the stage for increasing the grip, if the phone slides.
4. **Sample** – lay the prepared slide on the stage.
5. **Lighting** – Well align the light source with the Blips lens, under the sample. Use the *transparent foils* to lift the slide, for finding the right distance.
6. **Digitally zoom** the sample for a fine tuning; lock the automatic focusing.
7. Find tutorials, instructions, videos and our free **SMO App** at www.blips.pro